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Anleitung : **VE202601DE**

Operating Instructions LED Controller PLC300 – Easy

- 1 Safety instructions
- 2 Usage and functions
- 3 Connection options
- 4 Operation
- 4.1 Display & controls
- 4.2 Modes & functions
- 5 Troubleshooting
- 6 Environment and disposal
- 7 Warranty
- 8 Liability

IMPORTANT!
Read carefully before use! Store for later reference!

The content of this document has been checked for compliance with the products described. Nevertheless, deviations cannot be excluded, so that no guarantee can be given for full compliance. The manufacturer is not liable for damages due to missing or incorrect information. Changes or additions are always implemented without notice in the following issues.


1 Safety instructions
The instructions in all manuals must be followed. Failure to follow this instruction may result in property damage and personal injury or even mortal danger. With this controller, only the designated LED lights may be operated. (See → 2)
Switch off the LED controller before changing the pin assignment.

Never look directly into the LED bulbs when switching on or during operation!

2 Usage and functions
The controller is used only to connect and control Promicron LED spots and LED ring lights.

3 Connection Options
plug 3,5: A + B (see photo)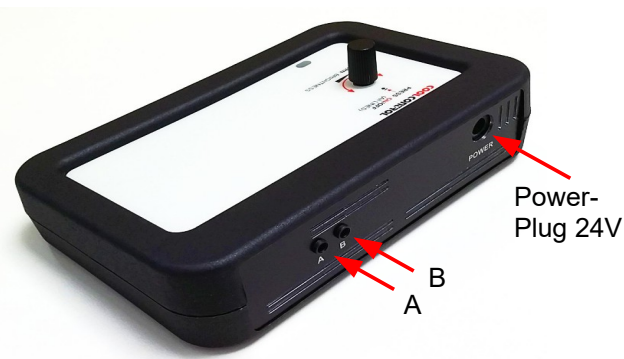

4 Operation

4.1 Display & controls
button / knob (see photo)

4.2 Modes & functions
Pushbutton / rotary knob **channel 1 (A+B)**
press : LED Controller ON / OFF
turn/rotate : brightness channel 1 (A+B)

5 Troubleshooting
Error : the LED lights are not working
Cause : spot defect or controller defective
Fix : please check spot on other channel; replace if necessary

6 Environment and disposal
The crossed-out wheeled bin symbol means that the product must be disposed of separately from household waste.

7 Warranty
We provide warranty within the scope of the legal provisions. Please return the device in the event of a fault together with a detailed description of the fault and return it to us with sufficient franking.

8 Liability
The product has been designed and manufactured for the intended use described in this document. Any other use is considered improper and may result in damage to the product or personal injury for which the manufacturer cannot be held responsible.

Manufacturer : Promicron AG, Bachmühlweg 24
74366 Kirchheim/Neckar, Germany

Instructions : **VE202601EN**